


# How To Crimp crucibles for Chip DSC

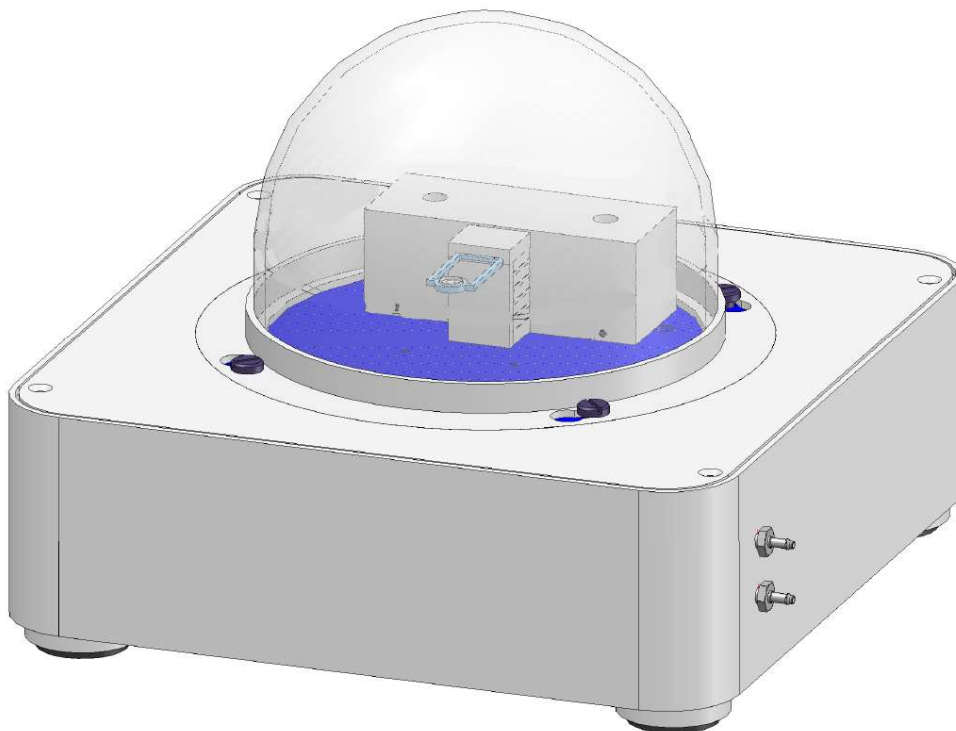

Linseis Messgeräte GmbH

Gerlach

Stand: 04.04.2019

## Index

1. The sample press and crimping tool.....	3
2. Prepare the Sample press and sample.....	4
4. Crimp the Crucible.....	5

**1. The sample press and crimping tool**




## 2. Prepare the sample press and sample

- Put the upper part from the crimping tool off



- Insert the crucible (make sure the bottom of the crucible is flat)

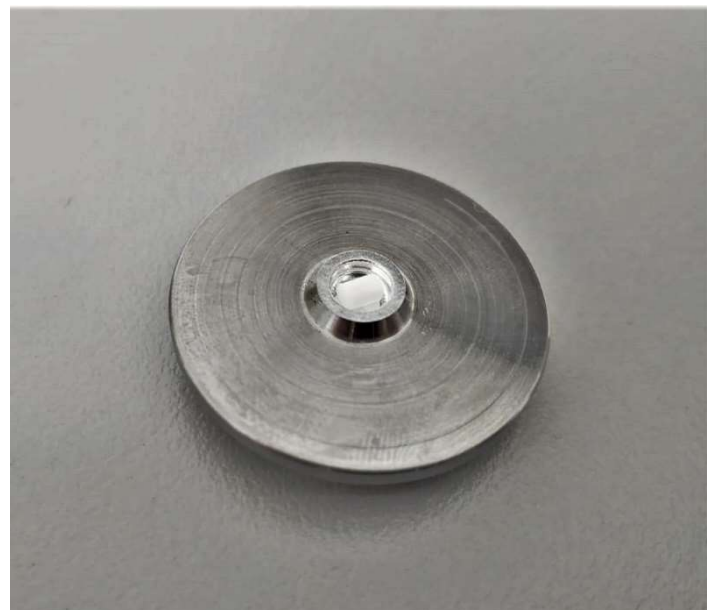

- Insert the sample into the crucible and put on the lid on the crucible

### 4. Crimp the Crucible

- Put the crucible holder back on the lower part of the crimping tool



- Pull down the rod of the crimping tool and crimp the crucible
- Use enough power on it, so that the crucible is pressed well



- Put the upper part of the crimping tool off from the sample press
- Your sample is ready now

